

Simultaneous Analysis of Quercitrin, Isoquercitrin, and Hyperoside on the Hitachi LaChrom Elite® Liquid Chromatography System

Hitachi High Technologies America, Inc.

Quercitrins modified with various carbohydrate glycosides via an ether linkage are often extracted from the hawthorn plant (*Hypericum japonicum*) for use as antioxidants¹. These flavinoid compounds, collectively known as quercitrins, might show health benefits in humans when dosed as supplements. Quercetin, itself, is a widely occurring compound, found in most produce, and the corresponding quercetin/saccharide ethers are more prevalent than free quercetin. Three common quercitrins—quercetin-3-O-galactoside (hyperoside, or HYP), quercetin-3-O-glucoside (isoquercitrin, or ISO), and quercetin-3-O-rhamnoside (quercitrin, or QUE)—were investigated for this application note. The goal of this study was to effectively resolve a mixture of the three components and show linearity and system suitability for each. Using the Hitachi LaChrom Elite® liquid chromatography system, these three compounds were easily resolved using a gradient pumping method and UV detection.


Experimental Conditions

Module	Conditions and Other
Pump (L-2130)	Mobile Phase A: 0.5% Acetic Acid Mobile Phase B: Acetonitrile See Gradient Profile Below
Autosampler (L-2200)	Injection Volume: 5, 3, and 1 μ L of a 16.7 μ g/mL solution of each component
Oven (L-2300)	30°C
UV Detector (L-2420)	UV λ : 355 nm
Column	Phenomenex® Luna ODS(2) 5 μ m 4.6X250 mm
Standards	QUE: Sigma-Aldrich® – Q3001, 90% pure ISO: BioChemika – 17793, >90% pure HYP: BioChemika – 83388, >97% pure Prepared at 50 μ g/mL each 50/50 MP A and B


Gradient Table for Quercitrins:

Time (min)	%A	%B	Flow Rate (mL/min)
0.0	84	16	1.5
15.0	80	20	1.5
15.1	84	16	1.5
20.0	84	16	1.5

Results - Linearity of Standards (HYP, ISO, and QUE, n=3):

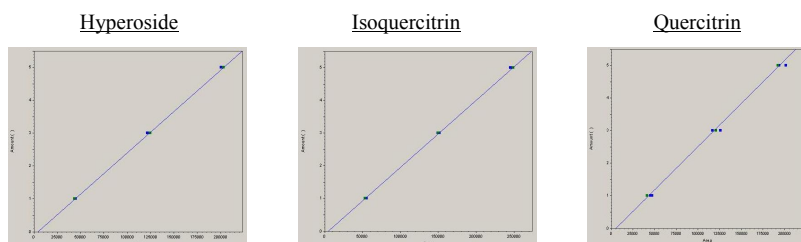

Results—Example Chromatogram of Quercitrins 3 μ L inj. Vol.

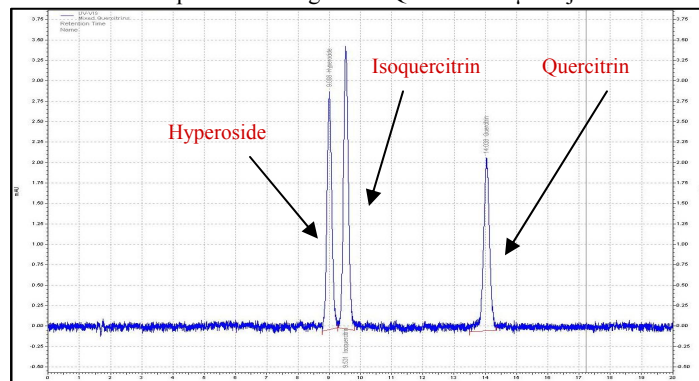

Reproducibility and Linearity of Standards

Compound	%RSD	Linearity
HYP 5 μ L inj.	0.6	$r^2 = 0.9999$
HYP 3 μ L inj.	1.2	-
HYP 1 μ L inj.	1.4	-
ISO 5 μ L inj.	0.7	$r^2 = 0.9999$
ISO 3 μ L inj.	0.6	-
ISO 1 μ L inj.	1.8	-
QUE 5 μ L inj.	2.5	$r^2 = 0.9991$
QUE 3 μ L inj.	2.8	-
QUE 1 μ L inj.	2.6	-

Discussion

Hitachi's LaChrom Elite® Liquid Chromatography System, equipped with a 5- μ m particle size Phenomenex® column, resolves the quercetin flavinoids hyperoside, isoquercitrin, and quercitrin. The system is suitable, and reproducibility of the standards is shown. Reproducibility (<3 %RSD) and system suitability (theoretical plates: $N > 1000$ and tailing factor: $T < 2.0$) are shown at the low quercitrins (HYP, ISO, and QUE) concentrations.

References:

1 – Chang, Q. et al. "HPLC Method for Simultaneous Determination of Hawthorn Active Components in Rat Plasma." *J Chromatogr B Biomed Sci Appl*, 688 (1) 1997.

*Hitachi is a registered trademark of Hitachi, Ltd. LaChrom Elite is a registered trademark of Hitachi High Technologies America, Inc. All other trademarks are the property of their respective owners.

Hitachi High Technologies America, Inc.

Life Sciences Division
 5100 Franklin Drive
 Pleasanton, CA 94588
 Toll Free: (800) 548-9001
 Email: Sales-LS@hitachi-hta.com
 Web site: www.hitachi-hta.com/hplc